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## Masson's Trichrome Stain Kit (Fast Green FCF Method)

**Cat:** G1343

**Size:** 7×50mL/7×100mL

**Storage:** RT, avoid light, valid for 1 year.

### Kit Components

Reagent		7×50mL	7×100mL	Storage
Reagent(A) :Weigert Hematoxylin Dyeing Solution	A1:Weigert Stain Solution A	25mL	50mL	RT, avoid light
	A2:Weigert Stain Solution B	25mL	50mL	RT, avoid light
Before use, mix with A1 and A2 in equal amount to form the Weigert Hematoxylin Dyeing Solution. It is not suitable to prepare it in advance.				
Reagent(B): Acid Ethanol Differentiation Solution		50mL	100mL	RT
Reagent(C): Masson Bluing Solution		50mL	100mL	RT
Reagent(D): Ponceau Fuchsin Solution		50mL	100mL	RT, avoid light
Reagent(E): Weak Acid Solution		50mL	100mL	RT
Reagent(F): Phosphomolybdic Acid Solution		50mL	100mL	RT, avoid light
Reagent(G): Fast Green Solution		50mL	100mL	RT, avoid light

### Introduction

In a narrow sense, connective tissue contains three kinds of fibers: collagen fibers, reticular fibers and elastic fibers. Collagen fibers are the most widely distributed and abundant fibers. Masson's Trichrome Staining is the most classical method in connective tissue staining. It is an authoritative and classical technique for collagen fibers staining. The Masson's Trichrome Staining usually refers to the staining of nucleus and the selective display of collagen and muscle fibers. The dyeing principle of this kit is related to the size of anionic dye molecules and the permeability of the tissues: the size of the molecules is reflected by the molecular weight, and the small molecular weight can easily penetrate the compact and low permeability tissues, while the large molecular weight can only enter the loose and high permeability tissues. However, the molecular weight of Fast Green is very large, so after Masson's Trichrome Staining, muscle fibers are red, collagen fibers are green (Fast Green), mainly used to distinguish collagen fibers from muscle fibers.

The characteristics of the Masson's Trichrome Stain Kit include: ◆ Stable staining; ◆ High fault tolerance, able to achieve satisfactory results within a large dyeing time span; ◆ Clear and bright colors; ◆ Widely used, suitable for staining tissues such as paraffin sections and frozen sections; ◆ The stained slices are stored for a long time and are not easily faded.

### Self Provided Materials

Fixative:10% formalin or 10% Neutral formalin., Series of ethanol, Distilled water.

### Protocol( for reference only)

#### For Paraffin Section

1. The recommended tissue thickness is 3-8μm. Dewax to distilled water.
2. (Optional)Before use, mix A1 with A2 in equal amount to prepare Weigert's Iron Hematoxylin Solution, stain with Weigert's Iron Hematoxylin Solution for 5-10mins.
3. (Optional)wash with distilled water to remove excess stain, differentiate with Acid Differentiation for 5-10s, wash with distilled water for 30s.
4. (Optional)Blue in Bluing Solution for 3-5mins. Rinse in distilled water for 30s.
5. Stain with Ponceau-Acid Fuchsin Solution for 5-10mins.
6. In the above step, mix distilled water with Weak Acid Solution in 2:1 ratio to prepare Weak Acid Working Solution. Rinse the section with Weak Acid Working Solution for 30s.
7. Discard the excess solution., differentiate in Phosphomolybdic Acid Solution for 1-2mins. Rinse the section with Weak Acid Working Solution for 30s.
8. Discard the excess solution, stain with Fast Green Solution for 1-2mins(See Note 11). Rinse the section with Weak Acid Working Solution for 30s.
9. Dehydrate quickly in 95% ethanol for 2-3s, then dehydrate with absolute ethanol twice for each time 5-10s.

10. Transparent in xylene for twice, each for 1-2min, seal with resinene.

#### For Frozen Section

1. The recommended tissue thickness is 8-12 $\mu$ m. After thawing and rehydrating the frozen sections, treat overnight with Bouin fixative at room temperature. (See Note 3)
2. Wash the slices with 75% ethanol until the yellow color completely fades away.
3. The subsequent operation is the same as steps 2-10 of paraffin section. (See Note 4)

#### Result

Collagen Fiber/Protein	Green
Cytoplasm, Muscle And Red Blood Cell	Red
Nucleus	Bluish Brown

#### Note

1. Slice dewaxing should be as clean as possible to avoid uneven color. The principle of this staining is related to the structure, and it is necessary to avoid non coloring caused by too thin tissue slices and overall darkening caused by too thick tissue slices. Therefore, it is recommended to have a slice thickness of 3-12 $\mu$ m.
2. The type of tissue and the choice of fixative have a significant impact on the staining results, so it is recommended to take a small number of sections for pre experiments for staining different samples.
3. If there is no Bouin fixative, G2310-Carnoy fixative II can be used instead for overnight treatment. After overnight treatment with Carnoy fixative II, soak twice in 75% ethanol and each for 1min, then conduct normal staining.
4. Frozen sections are more prone to dye adsorption due to the complete retention of cell components. Therefore, it is recommended to use Weak Acid Working Solution to dilute Fast Green Solution 2-4 times and stain for 0.5-1 minutes, followed by washing with Weak Acid Working Solution for 30s to avoid excessive blue staining.
5. This staining solution uses Weigert iron hematoxylin to stain the nucleus, which is a normal phenomenon that is inconsistent with the results of conventional hematoxylin staining. The purpose of staining is mainly to distinguish between collagen fibers and muscle fibers, and this staining step can generally be omitted.
6. The differentiation time of Acid Differentiation Solution should be determined according to the thickness of slice, the type of tissue and the old and new.
7. Masson Bluing Solution can also be replaced by G1865-Scott Bluing Solution or G1840/G1841-Lithium Carbonate Solution, 0.05%/1%.
8. Weak Acid Working Solution can make the color clearer and brighter. If it is used in a large amount, you can buy G2940-Weak Acid Solution or prepare 0.5-1% acetic acid solution to instead of Weak Acid Working Solution. If Fast Green is over dyed, you can replace the step "Rinse the section with Weak Acid Working Solution for 30s" by "Rinse the section with distilled water for 30s"
9. Phosphomolybdic Acid Solution is prone to failure and discoloration, and normally appears light yellow. If it has turned yellow green or blue during use, it is recommended to discard it.
10. Except for Reagent A and F, all reagents in this kit can be recycled for use, with a cycle count of 0.8-1ml for each section.
11. The time for Fast Green staining is adjusted according to the amount of collagen fibers in the tissue. Generally, tissues with more collagen fibers tend to stain faster, while tissues with less collagen fibers tend to stain slower and require longer staining time.
12. For your safety and health, please wear laboratory clothes and disposable gloves for operation.